# Bruksanvisning for 505CA-pumpehode

# 1 Innhold

---

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# EU DECLARATION OF INCORPORATION

1. Manufacturer: Watson Marlow Limited, Bickland Water Road, Falmouth, TR11 4RU, UK
2. This declaration of incorporation is issued under the sole responsibility of the manufacturer.
3. Object of the Declaration: 505CA.
4. The object of the declaration described above conforms in part with the relevant Union harmonisation legislation:  
*Machinery Directive 2006/42/EC*
5. The object of the declaration described above conforms in part with the following directive(s):  
*EMC Directive 2014/30/EU, RoHS Directive 2011/65/EU*
6. We undertake to transmit, in response to a reasoned request by the appropriate national authorities, relevant information on the partly completed equipment identified above. The method of transmission shall be by mail or email.
7. The pumphead is incomplete and must not be put into service until the machinery into which it is to be incorporated has been declared in conformity with the provisions of the Directive(s).

**SIGNED FOR ON BEHALF OF:**

Watson-Marlow Limited  
Falmouth, 31st December 2022

Simon Nicholson, Managing Director,  
Watson-Marlow Limited  
Watson-Marlow Fluid Technology Solutions  
Telephone: +44 (0) 1326 370370

A Spirax-Sarco Engineering plc company

**PERSON AUTHORIZED TO COMPILE THE TECHNICAL DOCUMENTS:**

Johan van den Heuvel

Managing Director

Watson Marlow Bredel B.V.  
Sluisstraat 7  
Delden  
Netherlands  
PO Box 47

Telephone: +31 74 377 0000

# UK DECLARATION OF INCORPORATION

1. Manufacturer: Watson Marlow Limited, Bickland Water Road, Falmouth, TR11 4RU, UK
2. This declaration of incorporation is issued under the sole responsibility of the manufacturer.
3. Object of the Declaration: 505CA.
4. The object of the declaration described above conforms in part with the relevant statutory requirements:

*Supply of Machinery (Safety) Regulations 2008, Electromagnetic Compatibility Regulations 2016, The Restriction of the Use of Certain Hazardous Substances in Electrical and Electronic Equipment Regulations 2012.*

5. We undertake to transmit, in response to a reasoned request by the appropriate national authorities, relevant information on the partly completed equipment identified above. The method of transmission shall be by mail or email.
6. The pumphead is incomplete and must not be put into service until the machinery into which it is to be incorporated has been declared in conformity with the provisions of the Directive(s).

**SIGNED FOR ON BEHALF OF:**

Watson-Marlow Limited  
Falmouth, 31st December 2022

Simon Nicholson, Managing Director, Watson-Marlow Limited  
Watson-Marlow Fluid Technology Solutions  
Telephone +44 (0) 1326 370370

A Spirax-Sarco Engineering plc company

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